

## Quick-guide for Pick&Mix microRNA PCR Panels

### Before starting the experiment

We recommend mixing your reagents with excess volume for pipetting. Typically, 10-25% excess is needed, depending on pipetting system. **Keep all reagents on ice while setting up.**

**Phase I: prepare RNA sample** (see Tips section page 51-52 for tips).

**Phase II (Steps 1-5): cDNA synthesis** (see page 45 for details).

- Adjust each template RNA to 5 ng/μl.
- For each sample, prepare and mix:

Reagent	Volume (μl) < 100 miRNA analyzed per sample	Volume (μl) > 100 miRNA analyzed per sample
5x Reaction buffer	2	4
Nuclease-free water	4.5	9
Enzyme mix	1	2
Synthetic spike in, optional replace with H <sub>2</sub> O if omitted	0.5	1
Template total RNA (5 ng/μl)	2	4
Total volume	10	20

Incubate for **60 min at 42°C** followed by heat-inactivation of the reverse transcriptase for 5 min at 95°C. Immediately cool to 4°C. If not used immediately, store at 4°C or freeze.

**Phase III (Steps 6-10): real-time PCR amplification** (see page 47 for details).

- Immediately before use, dilute your cDNA 100x. Suggestions for dilution procedure and volume needed for the specific Pick&Mix plate layout are given in the Protocol Table 10 (page 48). Add passive reference dye if recommended by instrument manufacturer.
- Mix 2x SYBR® Green MasterMix and diluted cDNA 1:1.
- Dispense 10 μl to each well of the 384-well PCR plate.
- Spin plate briefly in cooled centrifuge, wait 5 minutes while primers dissolve.
- Insert plate in cyclor, and run according to the following settings:

Polymerase Activation/Denaturation	95°C, 10 min
40 amplification cycles*	95°C, 10 s 60°C, 1 min, ramp-rate 1.6°C/s (100% standard on ABI instruments) Optical read
Melting curve analysis	

\*45 amplification cycles are required for LC480 instruments to allow collection of data for Cp values up to 40.

**Phase IV (Step 11): Data analysis**


ABI instruments: use manual baseline and threshold settings (see Tip 10, page 55).



## Workflow for Pick&Mix microRNA PCR Panels

### Phase I: Prepare RNA sample

See protocol page 51 for recommendations

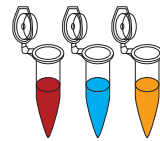

### Phase II: cDNA synthesis


See protocol page 45.

- Triplicate RT per sample is recommended
- a no enzyme RT negative control per study is recommended

### Phase III: real-time PCR amplification

See protocol page 47.

- Mix cDNAs with SYBR® Green
- Add cDNA:SYBR® Green mix to PCR plates

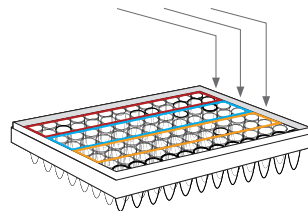


**ROX:** The SYBR® Green master mix, Universal RT does not include the ROX passive reference dye. Add as needed.

### Phase IV: Data analysis

Export data for further analysis:

- See data analysis guide online
- Data pre-processing, normalization and statistical analysis

**ABI 7900 and 7900 HT:** sds template files with pre-defined plate layout, cycling conditions and analysis settings are available at [www.exiqon.com/sds](http://www.exiqon.com/sds).

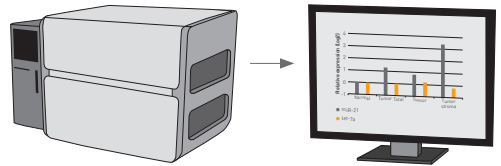

[exiqon.com/mirna-pcr](http://exiqon.com/mirna-pcr)

**EXIQON**  
Seek Find Verify

For life science research use only. Not for use in diagnostic procedures.

SYBR® Green is a registered trademark of Invitrogen. Concerning miRCURY LNA™ Universal RT microRNA PCR: NOTICE TO PURCHASER: LIMITED LICENSE Purchase of this product includes an immunity from suit under patents specified in the product insert to use only the amount purchased for the purchaser's own internal research. No other patent rights are conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.