

Quick-guide for individual microRNA LNA™ primer sets

Before starting the experiment

We recommend mixing your reagents with excess volume for pipetting. Typically, 10-25% excess is needed, depending on pipetting system. Prepare the **microRNA LNA™ PCR primer set**: Add 110 µl nuclease free water to each tube, vortex and spin down. Mix Fwd and Rev Primer 1:1 and spin down. Prepare the **Reference gene primer mix**: Add 220 µl nuclease free water, vortex and spin down. **Keep all reagents on ice while setting up.**

Phase I: prepare RNA sample (see Tips section, page 51-52 for tips).

Phase II (Steps 1-5): cDNA synthesis (see page 22 for details).

- Adjust each template RNA to 5 ng/µl
- Prepare and mix:

Reagent	Volume (µl)
5x Reaction buffer	4
Nuclease-free water	9
Enzyme mix	2
Synthetic spike in, optional replace with H ₂ O if omitted	1
Template total RNA (adjusted)	4
Total volume	20

Incubate for **60 min at 42°C** followed by heat-inactivation of the reverse transcriptase for 5 min at 95°C. Immediately cool to 4°C. If not used immediately, store at 4°C or freeze.

Phase III (Steps 6-10): real-time PCR amplification (see page 24 for details).

- When ready to set up the qPCR, dilute your cDNA 80x add 1580 µl H₂O pr 20 µl RT. Add passive reference dye if recommended by instrument manufacturer.
- To each well add:

Reagent	Volume (µl), 96/384 well plate, tubes or strips
SYBR® Green master mix	5
PCR primer mix	1
Diluted cDNA template	4
Total volume	10

Insert plate, tubes or strips in cyclor, and run according to the following settings:

Polymerase Activation/Denaturation	95°C, 10 min
40 amplification cycles*	95°C, 10 s
	60°C, 1 min, ramp-rate 1.6°C/s
	(100% standard on ABI instruments)
	Optical read
Melting curve analysis	

*45 amplification cycles are required for LC480 instruments to allow collection of data for Cp values up to 40.

Phase IV (Step 11): Data analysis

ABI instruments: use manual baseline and threshold settings (see Tip 10, page 55).



Workflow for individual primer sets

Phase I: Prepare RNA sample


See protocol page 51 for recommendations



Phase II: cDNA synthesis

See protocol page 22.


- Triplicate RT per sample is recommended
- a no enzyme RT negative control per study is recommended

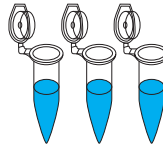

Phase III: real-time PCR amplification

See protocol page 24.

- Resuspend and mix PCR primers
- Add primers, cDNA:SYBR® Green mix to PCR plates

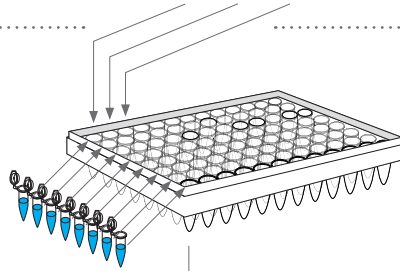


ROX: The SYBR® Green master mix, Universal RT does not include the ROX passive reference dye. Add as needed.

Phase IV: Data analysis

Export data for further analysis:

- See data analysis guide online
- Data pre-processing, normalization and statistical analysis

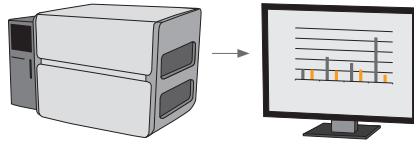

ABI 7900 and 7900 HT: 96- and 384-well sds template files with cycling conditions and analysis settings are available at www.exiqon.com/sds. Plate layout and detector must be added manually.

exiqon.com/mirna-pcr

EXIQON
Seek Find Verify

For life science research use only. Not for use in diagnostic procedures.

SYBR® Green is a registered trademark of Invitrogen. Concerning miRCURY LNA™ Universal RT microRNA PCR: NOTICE TO PURCHASER: LIMITED LICENSE Purchase of this product includes an immunity from suit under patents specified in the product insert to use only the amount purchased for the purchaser's own internal research. No other patent rights are conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.