

Guidelines for *in situ* hybridization on fresh frozen material using the miRCURY LNA™ ISH Optimization kits

These guidelines are a supplement to the miRCURY LNA™ microRNA ISH Optimization Kit (FFPE) manual V2.0. The manual can be downloaded from <http://www.exiqon.com/mirna-ish-kit>.

ISH protocols vary extensively due to different equipment set-up and Lab routines. This standard procedure guides the process of a manual *In Situ* Hybridization on human tissue using a DAKO Hybridizer (S2450).

Note that this protocol and the Exiqon miRCURY LNA™ microRNA ISH Optimization kits are not validated for use on Fresh Frozen material, and will likely require additional optimization. Please see page 23 -28 in the miRCURY LNA™ microRNA ISH Optimization Kit (FFPE) manual V2.0 for tips and troubleshooting.

General recommendations

Every step in the following procedure must take place in a clean and nuclease free environment. The operator must wear gloves during the entire process, and only use dry-sterilized glassware etc. (see below). in order to maintain RNase-free conditions. Use only ultrapure water or RNase free milli-Q H₂O in all reagents. All buffers/reagents applied in the assay must be RNase negative.

All glassware, including slides and coverslips, should be washed in a dishwasher. Wrap the items in aluminum foil so the surface or openings are covered and place a piece of indicator tape on the surface of the foil. Place items in an appropriate oven to be dry-sterilized for 8 hours at 180°C.

Note: for optimal *in situ* hybridization results on mouse tissue, it is recommended that the tissue specimens are obtained from perfusion fixed mice (see separate Guidelines for perfusion fixation). Before OCT-embedding, the fixed tissue samples must be cryo-protected with 20-30% sucrose in PBS overnight at 4°C. OCT-embedding/snap-freezing can be performed in isopentane at -80°C or liquid nitrogen. Store OCT-embedded samples at -80°C.)

microRNA ISH protocol

Before starting the experiment, please refer to the miRCURY LNA™ microRNA ISH Optimization Kit (FFPE) manual pages 12-18. Note the recommendations regarding optimization of the protocol using control probes. When using Fresh frozen samples, Step 1 (deparaffinization of FFPE slides) in the protocol is replaced by the following Steps:



Step 1a Cryo sectioning and fixation	Cut 10µm thick sections from OCT embedded FrFr blocks, mount them on a clean, non-contaminated, SuperFrost®Plus slides and let them air-dry for app. 15min (depending on the size of the tissue).Submerge the slides in 10% Neutral buffered Formalin (NBF) and leave it over night (o/n) at room temperature (RT).
Step 1b Wash	Transfer slides into container(s) with PBS and rinse for 3 x 3min (using fresh PBS for each wash).

Proceed to Step 2, page 19 of the miRCURY LNA™ microRNA ISH Optimization Kit (FFPE) manual V2.0.

Follow the rest of the protocol as described.

For life science use only. Not for use in diagnostic procedures.

Patents and Trademarks

Exiqon, AQ-Link, LNA, miRCURY, Hy3, Hy5, miRPlus, ProbeLibrary, ProbeFinder, and Genoview are registered trademarks of Exiqon, A/S, Vedbaek, Denmark.
Anthraquinone photochemistry products (AQ products) are covered by patents controlled by Exiqon A/S. Locked Nucleic Acids (LNA™) are covered by patents and patent applications owned by Exiqon A/S.

Disclaimers

Exiqon Products are for research use only and not for diagnostic or therapeutic use. The products may be used only for the buyer's internal research purposes and not for commercial use. The buyer may not resell products in their original or any modified form. The purchase of products does not include or carry an implied right or license for the buyer to use such products in the provision of services to third parties and a license must be obtained directly from Exiqon A/S for such use. DIG is licensed from Roche Diagnostics GmbH.

Exiqon A/S • Skelstedet 16 • 2950 Vedbaek • Denmark
Phone +45 45 650 929 • Fax +45 45 661 888
Exiqon, Inc • 12 Gill Street, Suite 1650 • Woburn, MA 01801 • USA
Phone +1 781 376 4150 • Toll free 888 647 2879
exiqon.com/contact • exiqon.com

