

In situ hybridisation on yeast with Locked Nucleic Acid in situ detection probes

Protocol prepared by Dr. Torben Heick, Department of Molecular Biology Aarhus University, Denmark

The protocol originates from the Singer laboratory (Long et al. (1997) Science 277, 383-387) and has been modified for yeast, *Saccharomyces cerevisiae*, cell cultures subjected to a heat-shock treatment. However, the outlined routines for fixation etc. should be applicable for yeast cell cultures in general.

Ref.: The protocol is modified from Thomsen et al. (2005) RNA, 11:1745-1748.

Fixation of yeast cells for RNA-FISH

- Cell cultures (10 mL) were grown at 25°C to an OD₆₀₀ of 0.1-0.3 and subjected to an instantaneous temperature-shift by addition of pre-heated media
- After incubation at the desired temperature and time, e.g. 31°C for 90 min, cells were fixed for 15 min at the experimental temperature by adding formaldehyde to a final concentration of 4%. This was followed by 30 min incubation at 20°C.
- Fixed cells were pelleted by centrifugation and washed three times in 1 ml wash buffer (1.2M sorbitol, 0.1M KH₂PO₄/K₂HPO₄, pH 6.5), and subsequently resuspended in 200-500 µL wash buffer.

Mounting on glass slides


- 10µl of the individual cell suspensions were plated on 14-well, 0.1% poly-L-lysine coated glass slides (Immuno-Cell Int.), washed in wash buffer containing 1% β-mercaptoethanol before spheroplasting for 10-15 min at 30°C in 10 µL oxalyticase solution (20mM vanadyl-ribonucleoside, 0.2% β-mercaptoethanol, 0.1 U/µL RNasin, 1.2M sorbitol, 0.1M KH₂PO₄/K₂HPO₄, pH 6.5, 0.1 mg/mL oxalyticase (Enzogenetics)).
- Subsequently, cells were washed for 5 min at 20°C; twice in wash buffer, once in (0.1M KH₂PO₄/K₂HPO₄, pH 6.5, 0.1% NP40) and once in (0.1M KH₂PO₄/K₂HPO₄, pH 6.5), and finally incubated with cold 70% ethanol for 15-30 min at -20°C.

Probe preparation

- Probe preparation was done by direct labelling of 10-20µg of 5'-end amino modified Locked Nucleic Acid in situ hybridization probe with 300µg of monofunctional Cy3 NHS-Ester (Amersham Pharmacia) in 0.1M NaHCO₃/Na₂CO₃, pH 9.0 overnight at 25°C in the dark. The probe was purified on a G-50 spin column and the concentration determined by OD measurement.
- For each well, 100ng of purified probe was mixed with 10µg salmon sperm DNA and 10µg yeast tRNA, lyophilized and resuspended in 5µL solution I (80% formamide, 10mM NaHPO₄ pH 7.0), denatured for 5 min at 95°C and finally mixed with 5µL of solution II (4xSSC, 20mM Vanadyl-ribonucleoside, 4µg/µL BSA, 0.1U/µl RNasin).

RNA-FISH

- Using a pipette, cells were drained for ethanol and washed in 10-20 µL (depending on well size) at 20°C, twice for 5 min in 2xSSC and once for 10 min in 40% formamide/2xSSC, 0.1% Triton X-100, before overnight incubation at 37°C with 10 µL of probe mix.
- Removal of probe solution was followed by washing steps in 10-20 µL volume: (i) twice in 40% formamide/2xSSC for 10 min at 37°C, (ii) once in 2xSSC/0.1% Triton X-100 for 10 min at 20°C, (iii) twice in 1xSSC for 10 min at 20°C, and (iv) twice in 1xPBS for 5 min at 20°C.
- Finally, 2.5µl of mounting solution (1xPBS pH 8.0, 80% glycerol, 0.1µg/mL DAPI) was applied to air-dried wells, which were subsequently covered with a cover slip and analysed by fluorescent microscopy.

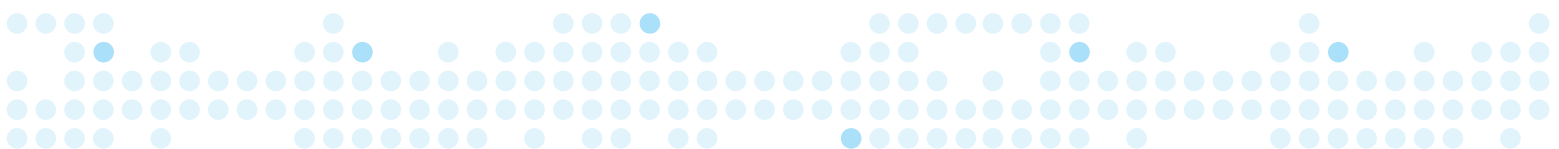

NB: Optimisation of the standard conditions by increasing hybridization stringency conditions can be achieved by increasing the formamide concentration (normally 40%) in hybridization- and all appropriate wash-buffers to 50%, 60% and 70%, respectively.

Imaging

Images were acquired on an Olympus BX51 microscope equipped with cooled Olympus DP50 CCD camera and analysis software.

Buffers

*Please note: For optimal fixation it may be critical to use fresh formaldehyde solutions. Fresh 4% solutions can be made from 16%, methanol free, formaldehyde or from solid paraformaldehyde (4% w/v).

For preparation of buffers please refer to :

Molecular cloning : a laboratory manual / Sambrook, Joseph; Russell, David W. --
3rd ed. -- New York: Cold Spring Harbor Laboratory, 2001.

"This protocol has been developed by a third party and not by Exiqon A/S or group companies ("EXIQON").

Thus, EXIQON cannot and will not warrant, represent or in any other way guarantee that the protocol and its content comply with your needs or expectations. EXIQON excludes all liability for any illegality arising from error, omission or inaccuracy in the protocol and takes no responsibility for the protocol or otherwise.

To the extent permitted by law, EXIQON excludes all liability in contract, tort (including negligence), breach of statutory duty or otherwise for any costs, losses, claims, damages, expenses or proceedings (including special, incidental or consequential loss or damage, loss of profits and wasted management time) incurred or suffered by you arising directly or indirectly in connection with the protocol and its content, including any loss, damage or expense arising from, but not limited to, any defect, error, imperfection, fault, mistake or inaccuracy with the protocol and its content. Any dispute relating to the protocol involving EXIQON shall be governed by Danish law."

