

# Quick-guide for Serum/plasma Focus microRNA PCR panels

(per panel, 2 x 96-well plates or ½ x 384-well plate)

## Before starting the experiment

Please note that this quick-guide refers to our Serum/plasma Instruction manual. We recommend mixing your reagents with excess volume for pipetting. Typically, 10-25% excess is needed, depending on pipetting system. Add this volume to the table volumes.

**Keep all reagents on ice while setting up.**

**Phase I: prepare RNA sample** (see Tips section, page 57-58 for tips).

**Phase II (Steps 1-5): cDNA synthesis** (see page 34 for details).

- For each sample, prepare and mix:

Reagent	Serum/plasma Focus Panel, Vol (µL)
5x Reaction buffer	4
Nuclease-free water	9
Enzyme mix	2
Synthetic RNA spike ins, optional replace with H2O if omitted	1
Template total RNA <sup>1)</sup>	4
Total volume	20

<sup>1)</sup> The equivalent of 16 µL original serum/plasma is used per 20 µL reverse transcription reaction.

Incubate for **60 min at 42°C** followed by heat-inactivation of the reverse transcriptase for 5 min at 95°C. Immediately cool to 4°C. If not used immediately, store at 4°C or freeze.

**Phase III (Steps 6-8): real-time PCR amplification** (see page 36 for details).

- Immediately before use, dilute your cDNA 50x: 980 µL water and 20 µL cDNA, and add 1000 µL 2x PCR Master mix. Add passive reference dye if recommended by instrument manufacturer.
- Dispense 10 µL to each well of the 96- or 384-well PCR plate(s) (two samples to a 384-well plate).
- Spin plate briefly in cooled centrifuge, wait 5 minutes while primers dissolve.
- Insert plate in cycler, and run according to the following settings:

Polymerase Activation/Denaturation	95°C, 10 min
40 amplification cycles*	95°C, 10 s 60°C, 1 min, ramp-rate 1.6°C/s (100% standard on ABI instruments) Optical read
Melting curve analysis	

\*45 amplification cycles are required for LC480 instruments to allow collection of data for Cp values up to 40.

## Phase IV (Step 9): Data analysis

ABI instruments: use manual baseline and threshold settings (see Tip 10, page 61).

# Workflow for Serum/plasma Focus microRNA PCR panels (per sample)

## Phase I: Prepare RNA sample

See protocol page 57 for recommendations



## Phase II: cDNA synthesis

See protocol page 34.

- Triplicate RT per sample is recommended
- a no enzyme RT negative control per study is recommended

*Illustration for a 96-well plate set-up*




## Phase III: real-time PCR amplification

See protocol page 36.

- Mix cDNAs with PCR Master mix
- Add cDNA:PCR Master mix to PCR plates

**ROX:** The ExiLENT SYBR® Green master mix, Universal RT does not include the ROX passive reference dye. Add as needed.

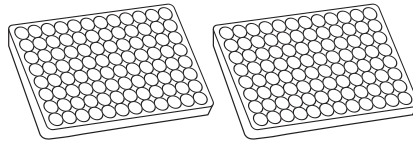


## Phase IV: Data analysis

Export data for further analysis:

- See data analysis guide online
- Data pre-processing, normalization and statistical analysis

**ABI 7500, 7500FAST, StepOne Plus, 7900 and 7900 HT:** 96- and 384-well sds template files with pre-defined cycling conditions and analysis settings are available at [www.exiqon.com/sds](http://www.exiqon.com/sds).

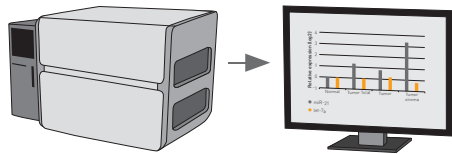

[exiqon.com/mirna-pcr](http://exiqon.com/mirna-pcr)

**EXIQON**  
Seek Find Verify

For life science research use only. Not for use in diagnostic procedures.

SYBR® Green is a registered trademark of Invitrogen. Concerning miRCURY LNA™ Universal RT microRNA PCR: NOTICE TO PURCHASER: LIMITED LICENSE Purchase of this product includes an immunity from suit under patents specified in the product insert to use only the amount purchased for the purchaser's own internal research. No other patent rights are conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.